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The screenshot displays the BioExpress dashboard interface. On the left is a dark blue navigation sidebar with the following items: Home, Devices, Alarms, Data, Tools, Setting, Account, and Technical Support. The main content area features a top navigation bar with a Home button, a notification area showing 'Device 2', 'SMS 50', and 'Voice 0', and a user profile for 'TEST01'. Below this is a breadcrumb trail for 'Home' and a yellow 'Notice' banner. The main section is titled 'Organization' and includes a 'Refresh' button and the text 'Update Time: 2023-12-26 07:49:26'. Two device status cards are shown: 'CryoMaster 6000_2' with a temperature of 9.3 °C and LN2 Level of 0 mm, and 'CryoMaster 6000_1' with a temperature of 8.4 °C and LN2 Level of 0 mm. Both devices are marked as 'Offline'.

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BioExpress

- Home
- Devices
- Alarms
- Alarm Project
- Alarm History**
- Data
- Tools
- Setting
- Account
- Technical Support

Home / Alarms / Alarm History

Device 1 SMS 25 Voice 0 test01

Home Devices Alarm Project Alarm History

Keyword: Enter Keyword Search Alarm Type: All Alarms Account Current Acco Alarm Time:

Select Alarm Time Range Search

Batch Processing Batch Delete

<input type="checkbox"/>	Device	Recipient	Alarm Time	Alarm Data	Operation
<input type="checkbox"/>	CryoMonitor 2000TL_1[2023120601]	CryoMonitor 2000TL-1_email,CryoMonitor 2000TL-1_SMS	2023-12-10 22:21:45	Temp:-182.9°C	Delete Processing
<input type="checkbox"/>	CryoMonitor 2000TL_1[2023120601]	CryoMonitor 2000TL-1_email,CryoMonitor 2000TL-1_SMS	2023-12-10 22:20:46	Temp:-182.9°C Liquid level:102ml	Delete Processing
<input type="checkbox"/>	CryoMonitor 2000TL_1[2023120601]	CryoMonitor 2000TL-1_email,CryoMonitor 2000TL-1_SMS	2023-12-10 22:19:46	Temp:-182.9°C Liquid level:102ml	Delete Processing
<input type="checkbox"/>	CryoMonitor 2000TL_1[2023120601]	CryoMonitor 2000TL-1_email,CryoMonitor 2000TL-1_SMS	2023-12-10 22:18:46	Temp:-182.9°C Liquid level:102ml	Delete Processing

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≡ Add User

i The Initial Password is 8888899999

* Username(Letters, Numbers, and Symbols @ .)

test01.2

* Name

test01.2

Phone

Please Enter the Phone Number

E-Mail

Please Input Your E-Mail

Select Organization

Select Organization

Select

Remark

Please Enter a Remark

≡ Menu Permission Settings [test01.1]

- System
 - Devices
 - Devices
 - Gateways
 - Alarm Devices
 - Alarms
 - Alarm Project
 - Alarm History
 - Datas
 - Daily Report
 - Search Data
 - Data Output
 - Data Chart
 - Tools
 - Sensor Tools
 - Gateway Tools
 - Alarm Tools
 - Setting
 - Login log
 - Operate log
 - Account
 - Change Password

≡ Operation Permission Settings [test01.1]

- System
 - Devices
 - Edit Device
 - View Data Chart
 - Print Data Chart
 - Manage Alarm Settings
 - treecode.caozuo.triggerdetailsedit
 - Alarms
 - Manage Alarm History
 - Delete Alarm History
 - Datas
 - Data Output
 - Search Data
 - Daily Report
 - Tools
 - Calibration Sensor

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